Safety Instructions for Puro[®] Printer YP-DP1

Safety instructions for Puro® Printer YP-DP 1

1 General safety information

NOTICE: Warning of damage to property and/or the environment.

The product was in perfect condition with regard to safety features when it left the factory. To maintain this condition and to ensure safe operation, the user must follow the instructions and observe the warnings in this manual.

Follow these safety precautions:

- Before operational startup, read the detailed documentation on <http://docs.puroscales.com/>.
- The equipment may only be used as intended for printing tasks.
- WARNING: Danger from incorrect supply voltage. It is essential that the power supply provided is used under all circumstances.
- CAUTION: Danger – print head may be hot! Allow the print head to cool down before replacing the paper roll.
- Observe the operating limits of the device.
- Position the device so that the power plug is freely accessible and the power cord does not present an obstacle or trip hazard.
- Unplug the power cord from the mains supply before cleaning.
- Make sure that no liquid enters the equipment.

2 Incoming goods inspection

Check the contents of the consignment for completeness. Check the contents visually to determine whether any damage has occurred during transport. If there are grounds for rejection of the goods, a claim must be filed with the carrier immediately. The sales organization must also be notified.

3 Before operational startup

NOTICE: Perform visual inspection.

Before operational startup as well as after storage or transport, inspect the product visually for signs of mechanical damage.

The product should not be put into operation if it displays signs of visible damage and/or is defective.

Access the detailed documentation by scanning the QR code

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Minebea Intec Bovenden GmbH & Co KG | Leinetal 2 | 37120 Bovenden